

## 10X Gmin Salts

- Begin with approx. 700 mLs MilliQ H<sub>2</sub>O
  - Weigh out and add ingredients with constant stirring:

▪ KH <sub>2</sub> PO <sub>4</sub>	47g
▪ K <sub>2</sub> HPO <sub>4</sub>	135g
▪ K <sub>2</sub> SO <sub>4</sub>	10g
▪ MgSO <sub>4</sub> -7H <sub>2</sub> O	1g
- Bring volume up to 1L
- Autoclave