

# YNB 4

## YNB (1x)

- Begin with approx. 800 mls of DI H<sub>2</sub>O
- Weigh out and add ingredients with constant stirring:
  - Yeast Nitrogen Base (without amino acids) 6.7g
- Bring up to 900 mLs with DI H<sub>2</sub>O
- Weigh out 20g of Glucose/Dextrose and add to 100 mLs of DI H<sub>2</sub>O
- Autoclave YNB and glucose separately
- Allow to cool in 50° water bath
- Combine YNB with glucose solution