

# YP 2

## YP (1.1x) (900 mL) Agar

- Begin with approx. 700 mls of DI H<sub>2</sub>O
- Weigh out and add ingredients with constant stirring:
  - Bacto Yeast Extract                    10g
  - Bacto Peptone                            20g
  
- Weigh out and add with stirring:
  - Bacto Agar                                20g
  
- Bring up to 900 mL with DI H<sub>2</sub>O
- Separate YP into two 1L bottles of 450 mL
- Autoclave
- Allow to cool in 50° water bath

*Verified 6/1/2005 by MaryAnn Hartman; Methods in Yeast Genomics, Cold Spring Harbor, 2000 ed.*