

S Buffer

- Begin with approx. 800 mLs of DI H₂O.
- Weigh out and add ingredients with constant stirring:
 - 1M KH₂PO₄ 43.55mL (final conc. of 43.55mM)
 - 1M K₂HPO₄ 6.45mL (final conc. of 6.45mM)
 - NaCl 5.85g
- Bring up to 1L with DI H₂O.
- Autoclave

Ian A. Hope (editor). *C. elegans: A Practical Approach*. Oxford University Press, New York, New York, 1999.